

## Condensation Particle Counter

Model 3786:RP-3786

Standard maintenance and calibration service for the Condensation Particle Counter Model 3786 consists of these procedures:

- 1) Review written instructions or explanation of problem from customer
- 2) Initial inspection
- 3) Update manual, firmware, and software as necessary
- 4) Clean saturator, growth tube, dryer, reservoir and optics
- 5) Inspect and replace as needed: wick, filters, orifices, screens, tubing, and o-rings
- 6) Flow Calibration
- 7) Check laser current and pump lifetime
- 8) Leak check
- 9) Zero check
- 10) Final electrical checkout and performance verification

Standard service includes minor repairs and replacement parts. Major repairs and replacement parts will be charged for parts and labor. TSI will contact the customer for approval before making any major repairs.

Turnaround time for standard service is approximately two weeks after receipt of the instrument. TSI service work is backed by a 90 day service warranty.

---

**TSI Incorporated** – 500 Cardigan Road, Shoreview, MN 55126 U.S.A

**USA** Tel: +1 800 874 2811

**UK** Tel: +44 149 4 459200

**France** Tel: +33 491 95 21 90

**Germany** Tel: +49 241 523030

**India** Tel: +91 80 41132470

**China** Tel: +86 10 8260 1595

E-mail: [particle@tsi.com](mailto:particle@tsi.com)

E-mail: [tsiuk@tsi.com](mailto:tsiuk@tsi.com)

E-mail: [tsifrance@tsi.com](mailto:tsifrance@tsi.com)

E-mail: [tsigmbh@tsi.com](mailto:tsigmbh@tsi.com)

E-mail: [tsi-india@tsi.com](mailto:tsi-india@tsi.com)

E-mail: [tsibeijing@tsi.com](mailto:tsibeijing@tsi.com)

Website: [www.tsi.com](http://www.tsi.com)

Website: [www.tsiinc.co.uk](http://www.tsiinc.co.uk)

Website: [www.tsiinc.fr](http://www.tsiinc.fr)

Website: [www.tsiinc.de](http://www.tsiinc.de)



TRUST. SCIENCE. INNOVATION.

Contact your local TSI Distributor or visit our website [www.tsi.com](http://www.tsi.com) for more detailed specifications.