

Survey of a Wake Field


Wake field measurements behind vehicles such as ships and aircraft, and objects such as airfoils and cylinders have been carried out to study energy dissipation, noise generation and suppression, interaction with propulsive devices and aerodynamic improvements. This type of information is generally obtained by mapping the flow field at various locations downstream of the vehicle or the body.

Fiberoptic LDV systems make this process of mapping the flow field by traversing the measuring volume a simple and straightforward process. The non-invasive nature combined with its ability to handle high turbulence and flow reversals makes the LDV technique ideally suited for these types of measurements.

The purpose of this experiment was to measure the wake field behind a two-dimensional airfoil. The flow field was surveyed to obtain velocity distribution and other flow properties in the wake region.

Experimental setup

A three component LDV system was used to measure the flow field behind the airfoil. The airfoil was mounted in a wind tunnel about a horizontal axis and set to have an angle of attack of about 5 degrees. The measurements were carried out with an argon-ion laser-based system using COLORBURST, COLORLINK, two 83 mm fiberoptic probes and IFA 755 Digital Burst Correlator. The measuring volume was traversed in a vertical plane to get the velocity variation.

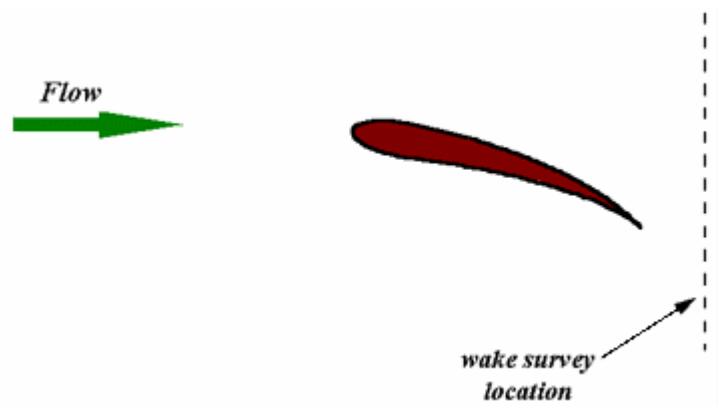


The free stream velocity of the air flow was about 9.5 m/sec. The flow was seeded using water droplets. The LDV system was operated in the coincidence mode to measure all three components of velocity simultaneously.

Results

The variation of the axial velocity mean (V_{mean}) plotted as a function of Z shows the velocity defect in the wake field behind the airfoil. All three components of velocity were also measured. The spanwise (w) and the vertical (u) components of velocity were significantly smaller than the streamwise velocity.

Mean Velocity & Re Stress Profile Flow behind an airfoil

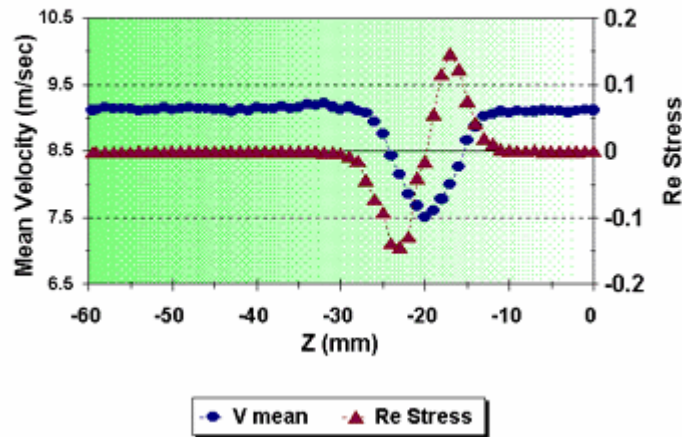

From these measurements all the higher order statistics, including Reynolds stress values, were calculated. The FIND-W software package was used for data acquisition and analysis. The figure also shows the variation of the Reynolds stress component ($u'v'$) contributed by the u' and the v' fluctuations.

Comments

These results show the ability of the LDV system to measure velocity very accurately so that mean velocities as well as higher order statistics (e.g., Reynolds stress) can be obtained accurately.



TSI Incorporated

Headquarters—Tel: +1 651 490 2811 **Toll Free:** 1 800 874 2811 **E-mail:** fluid@tsi.com

UK Tel: +44 1494 459200 **E-mail:** tsiuk@tsi.com **France** Tel: +33 491 95 21 90 **E-mail:** tsifrance@tsi.com

Germany Tel: +49 241 523030 **E-mail:** tsigmbh@tsi.com **Sweden** Tel: +46 8 595 13230 **E-mail:** tsiab@tsi.com

India Tel: +91 80 41132470 **E-mail:** tsi-india@tsi.com **China** Tel: +86 10 8260 1595 **E-mail:** tsibeijing@tsi.com

